# ***DrySyn*** ***Multi / Multi-M***

Betriebsanleitung  
Instruction Manual  
Mode d'Emploi  
Instrucciones de  
Empleo  
Istruzioni per l'uso

<b>D</b>	<b>DEUTSCH</b>	Seite	3 - 9
<b>E</b>	<b>ENGLISH</b>	page	10 - 16
<b>F</b>	<b>FRANCAISE</b>	page	17 - 23
<b>ES</b>	<b>ESPAÑOL</b>	página	24 - 30
<b>I</b>	<b>ITALIANO</b>	pagina	31 - 37

## INHALT

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## SUMMARY

STANDARD HARDWARE & OPTIONS.....	10
GENERAL INFORMATION .....	11
SAFETY INFORMATION .....	11
SETTING UP YOUR DRYSYN .....	12
SETTING UP YOUR REACTION .....	12
START YOUR REACTION HEATING.....	13
AFTER REACTION .....	13
CHOICE OF MAGNETIC STIRRING BARS / OPTIMUM SPEED.....	14
WARRANTY, LIABILITY & COPYRIGHT .....	14
QUESTIONS / REPAIR WORK.....	15

### STANDARD HARDWARE & OPTIONS

#### Standard Hardware

Product	Quantity	P/N
DrySyn MULTI	1	505-91900-00
DrySyn MULTI base plate	1	
DrySyn 25ml insert	3	
DrySyn 50ml insert	3	
DrySyn 100ml insert	3	
or		
Dry Syn MULTI-M		505-92600-00
DrySyn MULTI-M base plate	1	
DrySyn 100ml insert	3	
DrySyn 250ml insert	3	

## Options

Product	Quantity	P/N
Dry Syn MULTI-M 3-position Clamp	1	505-92090-00
Dry Syn MULTI 3-position Clamp	1	505-91090-00
Dry Syn MULTI 25ml Inserts	3	505-90020-00
Dry Syn MULTI 50ml Inserts	3	505-90030-00
Dry Syn MULTI 100ml Inserts	3	505-90040-00
Dry Syn MULTI-M 250ml Inserts	3	505-90050-00
Dry Syn MULTI reaction vial Inserts Sizes: 16,20 mm	3	505-90001-00
Dry Syn MULTI reaction vial Inserts, Sizes: 17,40 mm	3	505-90002-00
Dry Syn MULTI reaction vial Inserts, Sizes: 20,20 mm	3	505-90003-00
Dry Syn MULTI reaction vial Inserts, Sizes: 24,40 mm	3	505-90004-00
Dry Syn MULTI reaction vial Inserts, Sizes: 25,75 mm	3	505-90005-00
Magnetic stirring bars 15 x 6, (for 10ml), set of 3	1	509-53000-00
Magnetic stirring bars 25 x 10, (for 25 and 50 ml), set of 3	1	509-54000-00
Magnetic stirring bars 30 x 10, (for 100 and 250ml), set of 3	1	509-55000-00
Safety handle	1	505-91091-00

## GENERAL INFORMATION

-  **Unpack items carefully.**  
Inspect for damage and report such damage or missing parts to your supplier right away.
-  **Read your instruction manual carefully.** Take time to save time while working with your item. Make sure that every user has read and understood the instruction manual.
-  **Please store the instruction manual in a place easily accessible to every user.**

## SAFETY INFORMATION

-  **Please comply with all safety and accident-prevention regulations as in force for laboratory work !**
-  **Use extra care when working with flammable substances; refer to safety data sheets.**  
**Caution! Hazard of serious burns when running with hot plate over 50 °C.**
-  **Avoid touching the DrySyn or inserts at any time unless you are positive that it is cool.** The equipment will hold the heat for some time after it has been used.

## SETTING UP YOUR DRYSYN

4. Locate the DrySyn MULTI base plate on a hotplate stirrer (Maximum hotplate diameter 145mm).
5. Select the appropriate size of inserts for the flasks being used and carefully place them into the base plate, if you are only doing 1 reaction, use the central position for optimum stirring.
6. If available place the hotplate stirrer's temperature probe into the 3.5mm hole in one of the inserts.

## SETTING UP YOUR REACTION

### 1. INSPECT YOUR GLASSWARE

Please ensure your glassware is free of star cracks. Flasks can become significantly weakened around the perimeter by etching which occurs when the flasks are poorly loaded into glass washers. Please be aware that glassware with etching should not be used in the DrySyn.

### 2. ENSURE YOUR FLASK FITS PROPERLY

When selecting a flask, firstly ensure that the flask fits correctly by rolling it around in the insert. If the flask feels tight, or scratches the side of the DrySyn, or feels like it is sitting on an edge, the flask may be too big. Choose another flask. The DrySyn is manufactured to 0.2mm larger than the flasks supplied by Quickfit and Duran, please refer to the chart below.

DrySyn adapter size	Max flask diameter
25ml	41.2mm
50ml	51.4mm
100ml	65.2 mm
250ml	85.4 mm

3. Select a long thin stirring bar for optimum stirring. Carefully place the stirring bar into round bottomed flask. Do not drop stirring bars into the flask as this can cause breakage!

### 4. CLAMP YOUR FLASK SECURELY AT THE NECK

Although the flask is supported from beneath, it is also advised to clamp the neck to stop the flask from tilting. If using a condenser, it is essential to support the flask with a clamp to hold the weight of the condenser and reduce the pressure on the flask.

#### 5. USE A LAB JACK IF AVAILABLE

We strongly advise that you support the hotplate stirrer and DrySyn on a lab jack, this will offer the ability to lower the DrySyn away from the flask when cooling post synthesis as well as ensuring in an emergency the reaction flask can be lowered from the heat source. Lowering the DrySyn down by just 2-3mm during cooling will ensure that in the unlikely event that an oversize flask (See point 2) has been used it will not jam in the block.

#### 6. ENSURE THAT THE FLASK IS VERTICALLY POSITIONED IN THE INSERT

Flasks may not be spherical on the angle. You could increase stress on the flask if it is not vertically positioned.

7. Add a condenser or other glassware as desired. Solvents and reagents can be added more easily using a 2-neck flask.

8. Adjust the stirring speed to a suitable level for good mixing.

### START YOUR REACTION HEATING

1. Set the appropriate temperature on the temperature controller (if fitted) or via the hotplate control.

#### 2. DO NOT SET THE TEMPERATURE TOO HIGH

For low boiling solvents 5-10°C above the bpt is sufficient for reflux. For higher boiling solvents 10-20°C above the bpt will give good refluxing.

3. Ensure that there is adequate cooling water supply to the condenser to minimise loss of solvent.

#### 4. IF YOUR OPERATING TEMPERATURE IS ABOVE 150°C WE RECOMMENDED THAT YOU INSULATE THE FLASK WITH ALUMINIUM FOIL

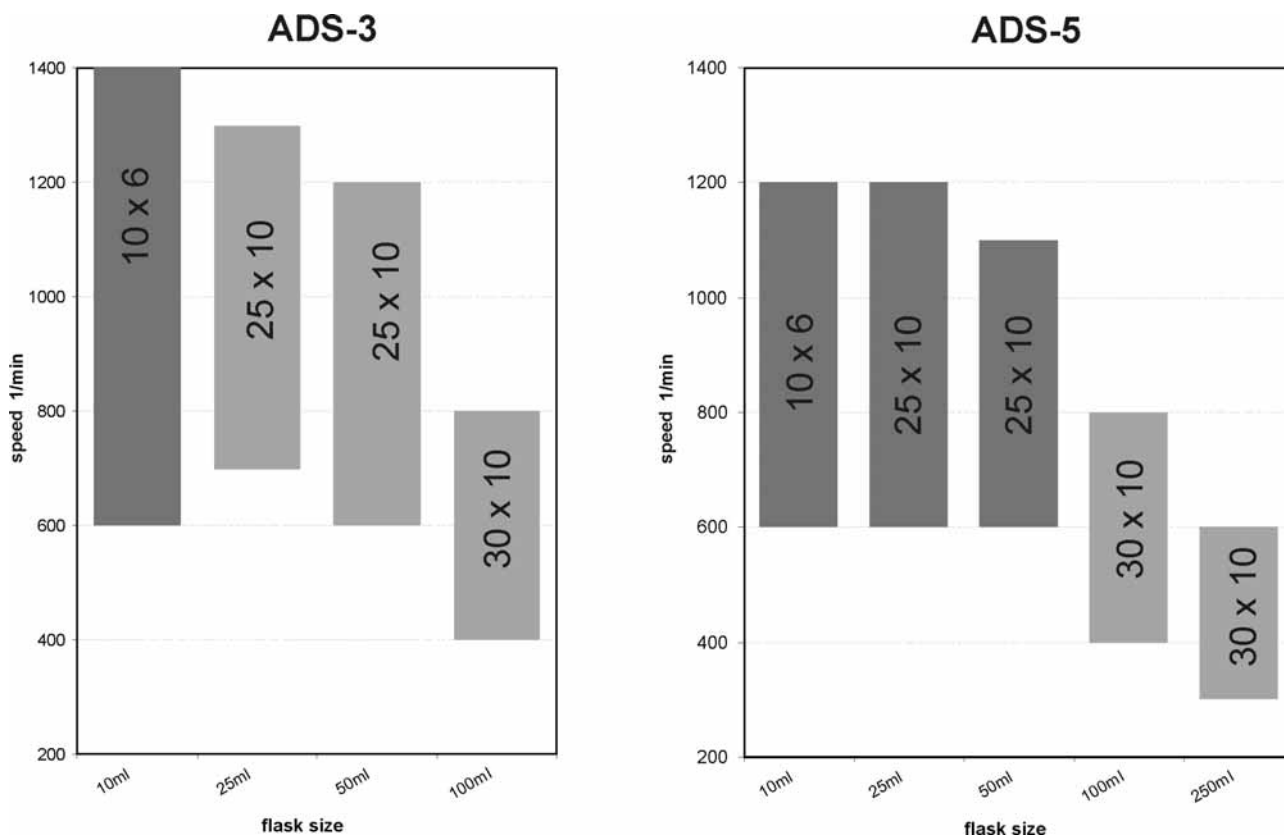
This reduces the thermal gradient across the glass and reduces stress on the flask at high temperatures. Important note! The maximum recommended temperature of the base for prolonged operation is 250°C. Exceeding this temperature may reduce the life of the stirring hotplate.

### AFTER REACTION

1. Please make your colleagues aware that the DrySyn could still be hot for some time after the hotplate has been switched off if you are leaving it unattended.

2. If available it is recommended that you lower the hotplate stirrer and DrySyn from the reaction flask on a lab jack during cooling.

## CHOICE OF MAGNETIC STIRRING BARS / OPTIMUM SPEED

Optimum speed settings of magnetic stirring bars (509-55000-00, 509-53000-00 and 509-54000-00, see chapter Standard Hardware & Options) measured in water.

## WARRANTY, LIABILITY & COPYRIGHT

### Warranty

Heidolph Instruments provides a three-year warranty on the products described here (with the exception of glass and consumable parts) if registered with enclosed warranty card or via internet ([www.heidolph.com](http://www.heidolph.com)). Warranty starts with the date of registration. Without registration warranty starts according to serial number. This warranty covers defects in materials and workmanship.

Transit damage is excluded from this warranty.

To obtain such warranty service, contact Heidolph Instruments (phone: +49 – 9122 - 9920-68) or your local Heidolph Instruments Dealer. If defects in material or workmanship are found, your item will be repaired or replaced at no charge.

Misuse, abuse, neglect or improper installation are not covered by this warranty promise.

Alterations to the present warranty promise need Heidolph Instruments' consent in writing.

### Exclusion Clause

Heidolph Instruments cannot be held liable for damage from improper use or misuse. Remedy for consequential damage is excluded.

### Copyright

Copyright in pictures and wording of the present instruction manual is held by Heidolph Instruments.

## QUESTIONS / REPAIR WORK

If any **aspect** of installation, operation or maintenance remains unanswered in the present manual, please contact us at the following address:

For equipment repairs jobs please call Heidolph Instruments (phone: +49 – 9122 - 9920-68) or your local authorized Heidolph Instruments dealer.

### Note

**You will receive approval for sending your defective item to the following address:**

Heidolph Instruments GmbH & Co. KG  
Lab Equipment Sales  
Walpersdorfer Str. 12  
D-91126 Schwabach / Germany  
Tel.: +49 – 9122 - 9920-68  
Fax: +49 – 9122 - 9920-65  
E-Mail: [sales@heidolph.de](mailto:sales@heidolph.de)

## SAFETY INFORMATION

When shipping items for repair that may have been contaminated by hazardous substances, please:

advise exact *substance*

take adequate *protective* action not to harm our parts receiving and service personnel

*mark* the pack IAW Hazardous Materials Act

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© Heidolph Instruments GmbH & Co. KG